

Hybridoma culture in the CellMaker: Practical production of monoclonal antibodies

Introduction

Antibodies are proteins within the immune system that identify and neutralise foreign proteins, or antigens. They exhibit a high level of specificity and affinity: two key functional characteristics that allow antibodies to be used as an integral tool in research, diagnostics, and therapeutics. Immunotherapy is gaining exceptional momentum in the scientific community, with a record number of antibodies being granted initial approval in the EU and US in 2018 (1). Compared to polyclonal antibodies, monoclonal antibodies (mAbs) tend to have better batch-to-batch consistency and less background reactivity, enabling them to be a useful tool for in vivo research and applications. According to a 2017 report, the monoclonal antibody market is projected to grow from \$403.7 million in 2017 to \$1,151.5 million in 2022 (2)— further exemplifying the magnitude and potential of this application. Hybridoma technology was the first mAb production method that allowed for in vivo scientific research and is still the primary and preferred platform for mAb derivatisation (3).



One of the most common and traditional ways to culture hybridoma cells utilises flasks stored in incubators. These can be labour intensive to prepare and maintain, and often yield inconsistent results due to hot and cold spots in incubators, high risk of contamination, and human error that comes with the workload of maintaining the culture.

Dr. Jean-Christophe Bourdon, Head of the p53 isoform laboratory at the University of Dundee, and his team have been using a CellMaker 8L system to effectively produce a range of p53 antibodies from murine hybridoma cells. In this experiment, batch cultures were run in both the CellMaker and flasks, and the respective production of two different mouse monoclonal antibodies were compared.

Methods

For this experiment, two production methods for two monoclonal antibodies were run in parallel: using a traditional flask method and using the CellMaker system. The method for each run is detailed below.

	Mouse Monoclonal Antibodies	
	C-alpha	mAb421
Seeded cell concentration	250,000 cells/ml	220,000 cells/ml
Flask volume	30ml	30ml
CellMaker volume	2L	2.3L
Duration of batch	7 days	4 days
Serum (FCS)	10%	10%
Media used	DMEM	DMEM

Flask Method

Under a fume hood, flasks were prepared and then stored in a 37°C incubator. They were checked routinely throughout each run to ensure no contamination was present.

CellMaker Method

The CellMaker 8L system that was used, enabled pH and dissolved oxygen monitoring and control. Under a fume hood, the sterile Cellexus bioreactor bag was fitted with two PTFE 0.2µm filters and an autoclaved pH probe. The empty bag was then fitted into the CellMaker enclosure, and the cells, media, and serum were pumped into the CellMaker bioreactor bag via a peristaltic pump and sterile tubing. For both runs, the following setpoints were defined by the user on the software:

Temperature	37°C
Airflow rate	0.1L per minute
pH	7.4
Dissolved Oxygen	99%

Temperature is controlled by an integrated Peltier heating system and monitored by a thermocouple in the enclosure. The pH is monitored via a Hamilton probe, and controlled by automatic addition of CO₂ to acidify the media when appropriate. A relative dissolved oxygen value is monitored by a fluorescence sensor built onto the bioreactor bag and controlled automatically by the software—introducing more air as necessary.


Downstream Methods

In this experiment, a protein G purification technique was used for both production methods. This can be substituted for peptide affinity chromatography or any other preferred purification technique.

Results

Results of total mAb produced are detailed in the table below. The amount of antibody produced from the total volume of the CellMaker was compared to the total volume that Dr. Bourdon would typically produce in one run with flasks: 10 flasks of 30ml each, equalling 300ml.

Monoclonal antibody	Production method	Total volume produced	Concentration mAb/ml	Total mAb produced
C-alpha	Flask	300ml	17.45µg/ml	5,235µg
	CellMaker	2000ml	13.41µg/ml	26,820µg
mAb421	Flask	300ml	16.67µg/ml	5,001µg
	CellMaker	2300ml	11.67µg/ml	26,841µg

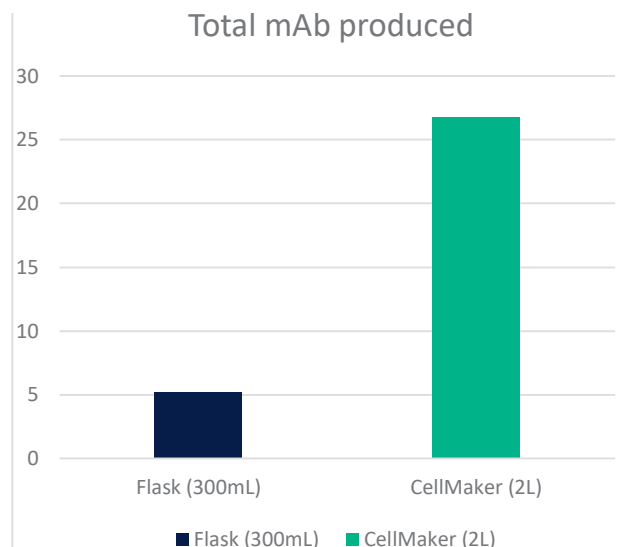

Discussion

The main advantages that the user observed using the CellMaker include the reduction in workload, small laboratory footprint, and stability. In terms of productivity and practicality, the CellMaker was the preferred method of the user compared to their traditional flask method.

Compared to other bioreactors, the CellMaker is unique in that it is the only single-use airlift bioreactor on the market. Airlift technology is desirable for mammalian culture due to the gentle yet efficient mixing technique. Conversely, powerful impellers in stirred tank bioreactors run the risk of shearing cell membranes. The disposable aspect of the CellMaker eliminates the need for sterilisation and validation between batches, while significantly reducing the risk of cross-contamination. The clear side panels of the CellMaker also allow for better visual checks than stainless steel systems.

Workload and space

The CellMaker comes in two sizes, the 8L and 50L. Using just the 8L system to its full working volume would equate to 267 flasks of 30ml. Such a large quantity of flasks would require significant workload and numerous hours to prepare and harvest. In addition, each of the 267 flasks should be checked throughout the week to ensure there is no contamination. This kind of workload and incubator storage space is not practical.

In contrast, the user noted that the CellMaker required approximately five minutes to prepare, three minutes to harvest, and no additional time in between to maintain. Checking for contamination was as simple as looking through the transparent bioreactor front panels. In addition, the 8L CellMaker can sit on a benchtop and does not require space in an incubator, as the Peltier element fitted to the enclosure is an accurate and precise way to control temperature.

Precise control and reproducibility

The CellMaker maintained constant temperature and pH throughout the run. Compared to incubator storage which can be contaminated easily and contain hot and cold spots, the contained, sterile CellMaker offers a much more stable environment for cells to grow and divide.

The media colour change due to lactic acid production is a visual indication that the run should be stopped, as antibody production is slowing down and the acidic environment runs the risk of degrading the quality of the antibody. Although this was observed in the flask cultures, the batch in the CellMaker exhibited different qualities. Throughout the runs, the pH was automatically maintained between values of 7.3 and 7.5, and the media did not change colour. This could indicate that the cells are remaining in normal metabolism, enabling long term growth and antibody production without increasing cost and labour. The resulting non-acidic environment could allow for greater quality of antibody.

The bag dynamically and tightly regulates oxygen, CO₂ and pH in the cell culture. Such parameters cannot be corrected in real-time in classic flask methods, preventing long term cell culture and limiting antibody production. The stability of the CellMaker enables long term cell culture and high titre antibody production.

Conclusions

For the same amount of time but far less workload and manual labour, the CellMaker provided a solution to scale up production of antibodies by five fold compared to flask production. Future experiments should test how long the culture could be sustained in the CellMaker before the media starts to acidify to evaluate the potential for even further antibody production. These results show that the CellMaker is a practical way to increase antibody production in a less labour-intensive manner in a single batch than the traditional method in multiple flasks.

“

We have found the CellMaker reliable, easy to use, low labour intensive and invaluable in our antibody production process. Our qualified team likes working with this bioreactor, using the advantages of disposable technologies and its precise regulation of biochemical parameters (pH, pO₂, temperature). At low-cost and with low labour, the CellMaker rapidly produces large quantities of high-quality antibodies with high specific activity. The CellMaker system is a vital part of the production process of our innovative and unique p53 antibodies.”

**Dr Jean-Christophe Bourdon, Principal investigator,
Head of p53 isoform research group, University of
Dundee**

References

1. Kaplon, H., & Reichert, J. M. (2018). Antibodies to watch in 2019. *MAbs*, 11(2), 219–238. doi: 10.1080/19420862.2018.1556465
2. Zaroff, S., & Tan, G. (2019). Hybridoma technology: the preferred method for monoclonal antibody generation for in vivo applications. *BioTechniques*, 67(3), 90–92. doi: 10.2144/btn-2019-0054
3. Kalia, R. (2017). Single Use Bioreactor Market Global Forecast to 2022. *Markets and Markets*, pp. 1–207.

Contact details

Cellexus International Limited
6 Riverside Court, Mayo Avenue
Dundee, DD2 1XD, UK

Email: enquiries@cellexus.com
Office: +44 (0) 1382 666357
www.cellexus.com

Contact us for more information or
to request a demonstration.